
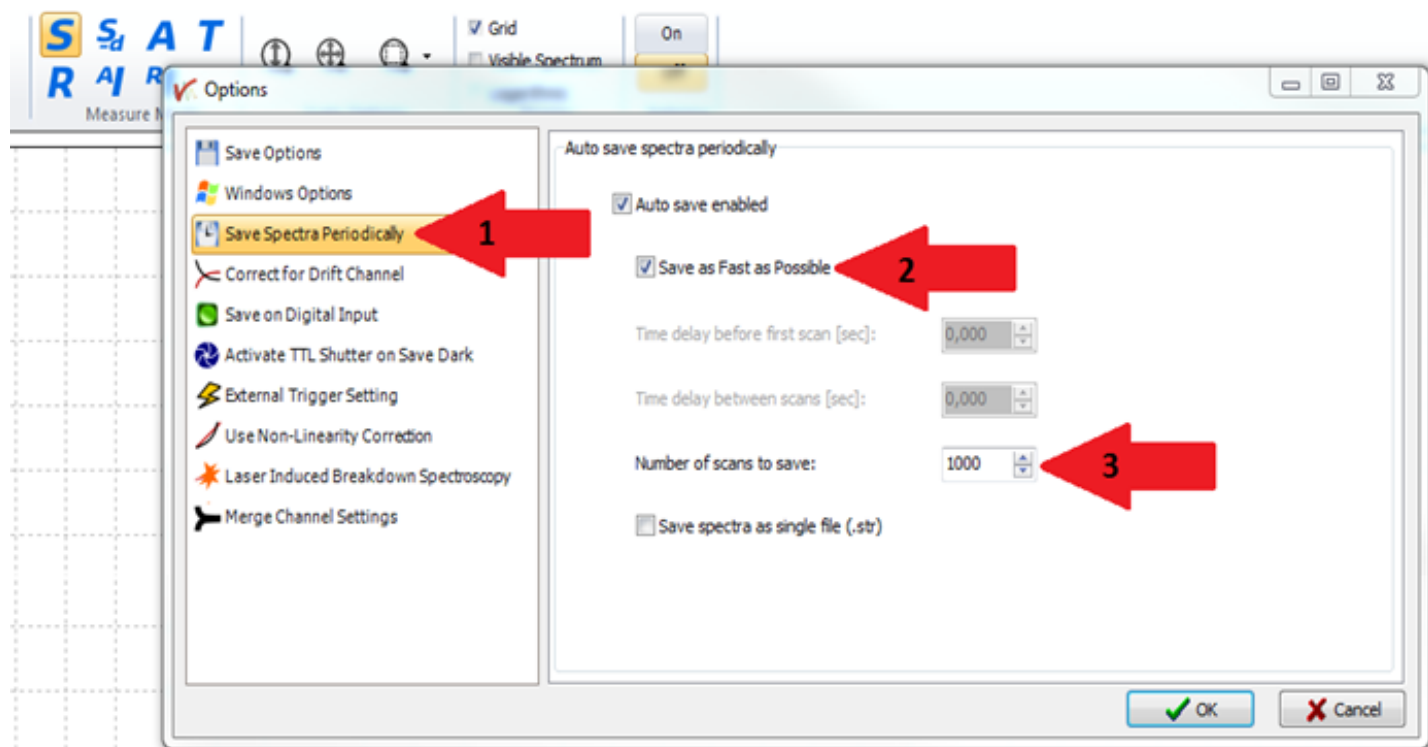



## How can I save my data periodically?

Please select Options from the toolbar. (See below picture).

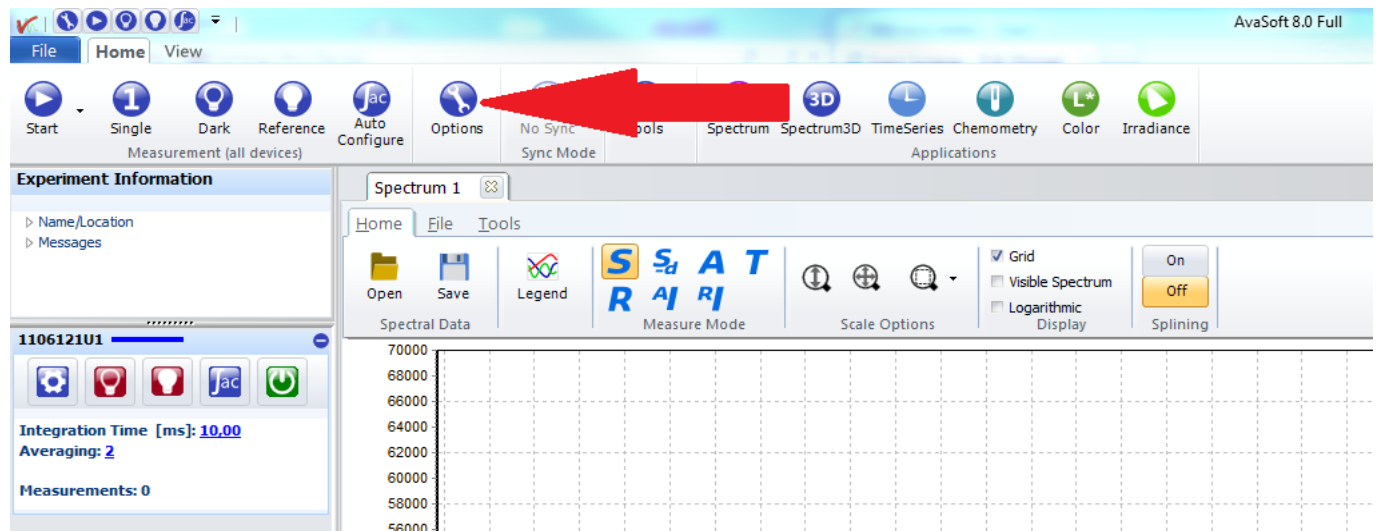

- 1) Please select "Save Spectra Periodically"
- 2) Please select from the option "Save as Fast as Possible".
- 3) Enter the number of scans you need to save.

After your measurements your data has been saved in different AvaSoft 8 files.

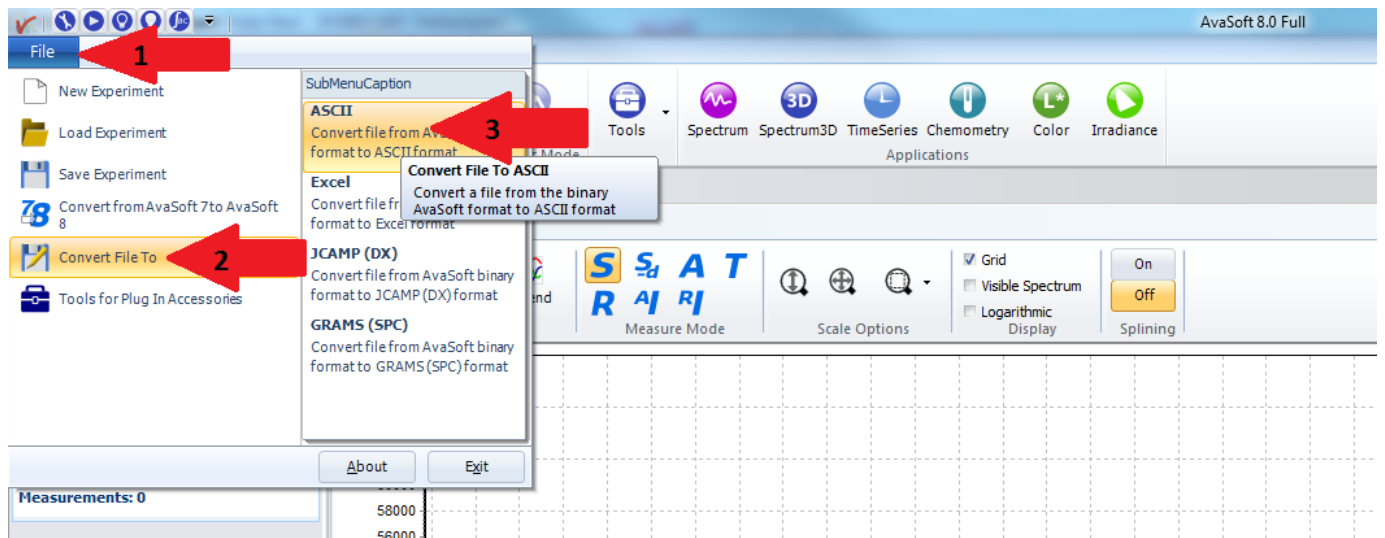

To convert these files to ASCII files you need to do the following.

- 1) Go to "File"
- 2) Go to "Convert file to"
- 3) Select "ASCII"

Select the files you need to convert to ASCII files.

Now your files have been converted to ASCII files.